


Seg-Flow® 4800 Tubing Replacement Schedule

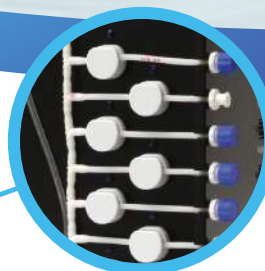

Seg-Flow Manifold
1, 2




Seg-Flow Pump*
3a, 4a

A little care goes a long way.

To keep your Seg-Flow® 4800 system at its best, Flownamics recommends periodic maintenance of the system. This will help to avoid issues such as tubing blockage, fluid delivery failure and incorrect fluid sensor readings that could prevent the Seg-Flow from successfully delivering a sample during a run.



5 Vessel Tubing Assemblies

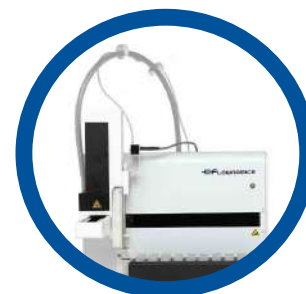

FlowFraction 400 7



9 Sample-Mod 300



6 Reagent/Waste Bottle



FlowFraction 400 Wash Pump 8

For quantities and pricing, contact your local distributor or manufacturer's rep. Or contact us directly at 800.932.6989.

Part Numbers & Schedule - Flownamics Products

	Part Number	Description	Replacement
1	100-1168	Seg-Flow® 4800 manifold tubing (without Reagent-Mod)	Every 3 months
2	100-1170	Seg-Flow 4800 manifold tubing (with Reagent-Mod)	Every 3 months
3a	100-1113	Seg-Flow 4800 sample delivery pump tubing	Every 3 months
4a	100-1114	Seg-Flow 4800 waste delivery pump tubing	Every 3 months
5	100-1143 100-1153	Vessel tubing assemblies (all variations)	Every bioreactor run
6	100-1098	Reagent and waste bottle tubing	Immediately if there is visible blockage/damage. If not, replace with annual PM.
7	100-1099	FlowFraction™ 400 tubing	Every 3 months
8	100-1111	FlowFraction 400 wash pump tubing	Every 6 months
9	100-1109	Sample-Mod 300 tubing	Every 3 months
*If your Seg-Flow 4800 has the first generation cassette pump, the following part numbers replace 3a and 4a from above.			
3b	100-1092	Seg-Flow 4800 sample delivery pump tubing (cassette pump)	Every 3 months
4b	100-1093	Seg-Flow 4800 waste delivery pump tubing (cassette pump)	Every 3 months



Part Numbers & Schedule - Integrations

	Part Number	Description	Replacement
10	100-1101	BioProfile® FLEX interface tubing	Every 3 months
11	100-1102	Cedex HiRes interface tubing	Every 3 months
12	100-1106	Cedex Bio HT interface tubing	Every 3 months
13	100-1108	BioProfile FLEX2 interface tubing	Every 3 months
14	100-1110	Kaiser Raman Rxn2™ interface tubing	Every 3 months
15	100-1119	BioProfile FLEX2 and FlowFraction tubing without Sample-Mod	Every 3 months
16	100-1094	YSI interface tubing	Every 3 months
17	100-1095	BioProfile 100, 300, 400 interface tubing	Every 3 months
18	100-1096	Vi-CELL® XR interface tubing	Every 3 months
19	100-1118	Seg-Flow pump tubing for BioProfile 100, 300, or 400	Every 3 months

Flownamics, Inc.

800-932-6989

contact@flownamics.com | www.flownamics.com